



Link2 LIVE Rad Trolley from Phoenix Dosimetry



The trolley comes ready assembled.

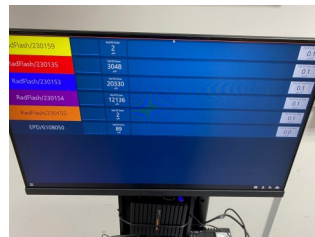


Step 4: Make sure the monitor is switched on and then wait for the software to boot up and the dose / doserates are displayed alongside the colour coded EPD's.

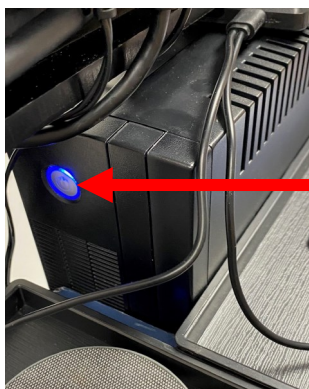


Step 1: Connect the 13 amp power cord to a 240v supply and then power on the UPS via the blue button, make sure the Radflash EPDs have the Bluetooth enabled by holding down the button on the back until it flashes blue.



Step 5: To reset the dose hold the EPD very close to the link 2 and this screen appears. Then just follow the instructions keeping the Radflash close.

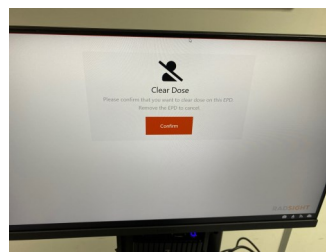

Step 2: Place the Radflash onto the charging stations, when charging they will go RED. Make sure they are dead centre or they will not charge—See red one not charging. Note: sometimes the clips interfere with the charging - if this is the case please remove the clip.

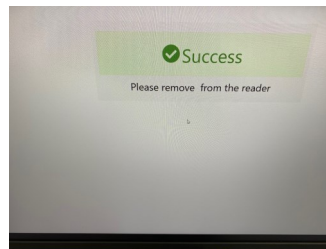

Step 6: Confirm with the mouse and the cumulative dose is now reset and the screen will default back.



Step 3: Switch on the Link2 via the blue button.